## BC-521 Gadolinium Loaded Liquid Scintillators

BC-521 is formulated to yield the highest light output possible as well as long term chemical stability. The standard gadolinium concentration is 0.5% by weight, but other concentrations up to 1.5% can be supplied. Since the liquid is normally used in large tanks containing several hundred liters, it employs a high flash point solvent for safety purposes.

The principal applications of BC-521 are for neutron spectrometry and neutrino research. It is ideal for use in large tanks for neutron multiplicity experiments.

Gadolinium has the highest thermal neutron capture cross-section of any element. The neutron capture reaction yields beta particles and several gamma rays having a total energy of about 8 MeV. Delayed coincidence and pulse shape discrimination techniques can be employed.

Scintillation Properties	
Light Output, %Anthracene	68
(for 1% Gd concentration), % Anthracene	57
Wavelength of Maximum Emission, nm	424
Decay Time, short component, ns	3.6
Bulk Light Attenuation, meters	>4
Atomic Composition	
No. of H Atoms per cc (x10 <sup>22</sup> )	5.25
No. of C Atoms per cc (x10 <sup>22</sup> )	4.00
No. of Electrons per cc (x10 <sup>23</sup> )	2.97
Ratio H:C Atoms	1.314

## General Technical Data -

Density	0.89 g/cc
Refractive index	1.50
Flash Point	44°C
Gadolinium Content	0.5%, w/w



## BC-521 Gadolinium Loaded Liquid Scintillator



## **Emission Spectrum**



Saint-Gobain Crystals

www.crystals.saint-gobain.com

Manufacturer reserves the right to alter specifications. ©2005-2018 Saint-Gobain Ceramics & Plastics, Inc. All rights reserved.